# Protocol for Real-Time RT-PCR assay with Lyoph P&P CHIKV Duo (adapted to TaqMan platform)

### For RUO (Research Use Only)

#### Vial content

Lyoph P&P CHIKV Duo contains lyophilized primers and probe (P&P) for detection of chikungunya virus (CHIKV). It must be used in combination with a generic RT-qPCR kit.

#### Design

RT-qPCR systems included in Lyoph P&P CHIKV Duo:

- Development of a TaqMan RT-PCR assay without RNA extraction step for the detection and quantification of African Chikungunya viruses. Pastorino et al., J Virol Methods, 2005. <u>DOI</u>: 10.1016/j.jviromet.2004.11.002
- Chikungunya fever in travelers returning to Europe from the Indian Ocean region, 2006. Panning et al., Emerg Infect Dis, 2008. <u>DOI</u>: 10.3201/eid1403.070906

#### Caution

Vials containing primers and probe mix (Lyoph P&P) must be stored at -20°C in the dark after reception. Stable for at least 6 months under the described conditions.

#### Instructions

This protocol is adapted to the following kit and RT-qPCR platform:

- SuperScript™ III Platinum™ One-Step qRT-PCR Kit (ThermoFisher)
- CFX (BIORAD) RT-qPCR thermal cycler

Use of other kits or RT-qPCR platforms may require adaptation of the protocol.

### I. Rehydration of Lyoph-P&P

- Write the date on the vial before opening
- Resuspend Lyoph P&P as follows (these steps are critical to ensure adequate homogenization):
  - ✓ Add nuclease-free water or RE buffer as described in Table 1
  - ✓ Homogenize by pipetting 10 to 20 times the volume up and down in the glass vial
  - ✓ Incubate rehydrated P&P at room temperature for 10 min
  - ✓ Perform a second series of 10-times pipetting.

**Table 1.** Lyoph-P&P regeneration; the number of tests/vial is indicated on the Lyoph P&P vial label.

Packaging (number of tests/vial)	8+1	16+2	24+2	32+3	48+3	96+5
H₂O or RE buffer (μL)	63	126	182	245	357	707

## II. <u>Preparation of the PCR reaction mix (for SuperScript™ III Platinum™ One-Step qRT-PCR Kit - ThermoFisher)</u>

 Table 2. Preparation of the PCR mix. \*Rehydration performed as indicated in Table 1.

	Volume/PCR test (μL)
2X Reaction mix	12.5
Rehydrated Lyoph P&P*	7
SSIII/Taq enzyme mix	0.5
Total (1 PCR test)	20
Template RNA (H₂O for Ctrl-)	5
Final volume	25

# III. Cycling program and RT-qPCR (for SuperScript™ III Platinum™ One-Step qRT-PCR Kit - ThermoFisher)

50°C	15min		
95°C	2min		
95°C	15sec	45 cycles	
60°C	30sec (plate read)	45 Cycles	

Read: FAM-QSY

Rehydrated Lyoph-P&P can be stored at 4 °C in the dark for up to 7 days.